



**Kalrez<sup>®</sup>** perfluoroelastomer  
parts

From DuPont Performance Elastomers

## Kalrez<sup>®</sup> perfluoroelastomer seals developed specifically for couplings according to DIN11864-2 Form A

Couplings based on DIN11864-2, Form A, assume complete filling of the groove after compression of the o-ring in order to form a hygienic seal (Figure 1). Exposure of the coupling to various chemicals, cleaning agents, steam and elevated temperatures, can cause the o-ring material to swell, resulting in intrusion of the o-ring into the process stream (Figure 2). Consequently, a dead space may be formed, leading to contamination and loss of the hygienic properties of the coupling. Additionally, the o-ring may be damaged, break off and contribute to particulates in the process.

Some elastomeric compounds, like EPDM and fluoroelastomers, have limited chemical and thermal resistance, especially when exposed to the CIP and SIP cycles being used in today's pharmaceutical and biopharmaceutical processes. Kalrez<sup>®</sup> perfluoroelastomer parts, with their superior chemical resistance, provide materials that can avoid these potential volume swell problems and are being used more often in today's environments.

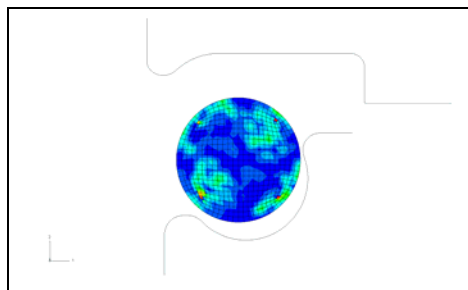


fig. 1 : O-ring section before installation

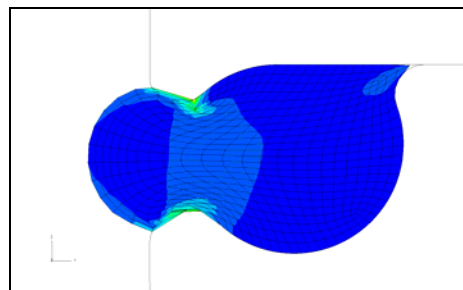


fig. 2 : intrusion into the process stream

Kalrez<sup>®</sup> parts, however, do exhibit thermal expansion - a physical property that cannot be avoided and could also lead to intrusion in the process stream. Engineers at DuPont Performance Elastomers S.A in Geneva created a new design with the help of Finite Element Analysis (FEA).

Our new design is a D-shape (Figure 3) and precisely sized to create sufficient free volume in the groove. This allows the Kalrez<sup>®</sup> material to expand without intrusion into the process stream (Figure 4).

<sup>1</sup>Kalrez<sup>®</sup> parts based on compounds 6230 (black) and 6221 (white) are well suited for these types of designs and couplings and comply with the requirements of U.S. FDA regulation 21 CFR 177.2600 and the extractive requirements of 21 CFR 177.2400(d)(1). They have also been tested and in accordance with requirements of a United States Pharmacopeia Class VI testing protocol and meet the requirements of USP Class VI polymer.

<sup>1</sup>Kalrez<sup>®</sup> perfluoroelastomer parts are not routinely tested using the USP testing protocol. Cured samples made only from compounds 6221 and 6230 have been tested in accordance with USP protocols and meet the requirements of a USP Class VI polymer. USP testing was done to support use of Kalrez<sup>®</sup> parts in pharmaceutical processing and food processing applications. While USP Class VI compliant materials are not required for pharmaceutical and food processing applications, many pharmaceutical and food processing customers, including customers seeking ISO9000 certification, have requested compliance. Testing of any finished article that incorporates Kalrez<sup>®</sup> perfluoroelastomer parts is the responsibility of the manufacturer or seller of the finished article if certification that meets USP standards is required.

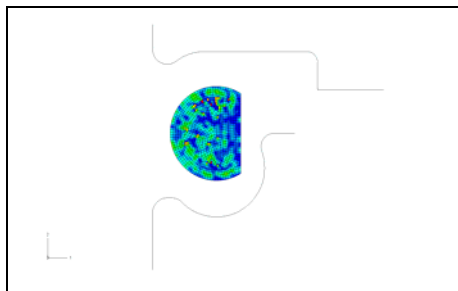


fig. 3 : Kalrez® D-shaped section ring

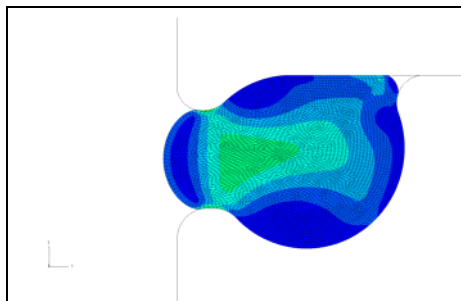


fig 4. : Kalrez® D-ring does not intrude in the process stream

A major French pharmaceutical company has successfully tested our design. The photo in Figure 5 shows a typical material that has swollen and taken a permanent deformation after only 10 SIP cycles – the intrusion into the process stream is obvious. Figure 6 shows the Kalrez® design after 22 SIP cycles, where swelling is minimal along with minimal intrusion into the process stream.



Kalrez® perfluoroelastomer seals for DIN11864-2 are available in nominal dimensions DN10, DN15 and DN20. Other dimensions can be made available upon demand.

For further information please contact one of the offices below, or visit our website at [www.dupontelastomers.com/kalrez](http://www.dupontelastomers.com/kalrez)

**Global Headquarters – Wilmington, DE USA**

Tel. +1-800-853-5515  
+1-302-792-4000  
Fax +1-302-792-4450

**European Headquarters - Geneva**

Tel. +41-22-717-4000  
Fax +41-22-717-4001

**South & Central America Headquarters - Brazil**

Tel. +55-11-4166-8978  
Fax +55-11-4166-8989

**Asia Pacific Headquarters - Singapore**

Tel. +65-6275-9383  
Fax +65-6275-9395

**Japan Headquarters – Tokyo**

Tel. +81-3-6402-6300  
Fax. +81-3-6402-6301

The information set forth herein is furnished free of charge and is based on technical data that DuPont Performance Elastomers believes to be reliable. It is intended for use by persons having technical skill, at their own discretion and risk. Handling precaution information is given with the understanding that those using it will satisfy themselves that their particular conditions of use present no health or safety hazards. Since conditions of product use and disposal are outside our control, we make no warranties, express or implied, and assume no liability in connection with any use of this information. As with any material, evaluation of any compound under end-use conditions prior to specification is essential. Nothing herein is to be taken as a license to operate or a recommendation to infringe on patents. While the information presented here is accurate at the time of publication, specifications can change. Check [www.dupontelastomers.com](http://www.dupontelastomers.com) for the most up-to-date information.

**Caution:** Do not use in medical applications involving permanent implantation in the human body. For other medical applications, discuss with your DuPont Performance Elastomers customer service representative and read Medical Caution Statement H-69237.

DuPont™ is a trademark of DuPont and its affiliates.

Kalrez® is a registered trademark of DuPont Performance Elastomers.

Copyright © 2005 DuPont Performance Elastomers. All Rights Reserved.